



PREPOR GP Filter Cartridges

- liquid filters
- glass microfibre / polypropylene

PREPOR GP is a new prefilter that combines the strength of polypropylene with the microbial retention of glass fibre for demanding applications such as long term exposure to steam, high differential pressures or aggressive chemicals.

The combined media will also provide a significant microbial reduction that makes PREPOR GP equally suitable for bioburden reductions in pharmaceutical liquids as well as offering excellent protection to sterilizing grade membrane cartridges. By using graded density media, PREPOR GP has a higher voids volume (95%) and greater dirt holding capacity than surface filtration membranes which means that filtration costs are reduced without affecting the product quality. PREPOR GP can also provide excellent prefiltration to membrane filters in proteinaceous and high contamination applications by extending the life of the membrane cartridge and hence reducing filtration costs.


Features and Benefits

- Combined media for microbial retention and mechanical strength
- Graded density media gives increased dirt holding capacity
- Suitable for bioburden reduction and fine prefiltration
- Pleated construction with rigid core and sleeve




Note: PREPOR is a registered trademark of Parker domnick hunter

Performance Characteristics

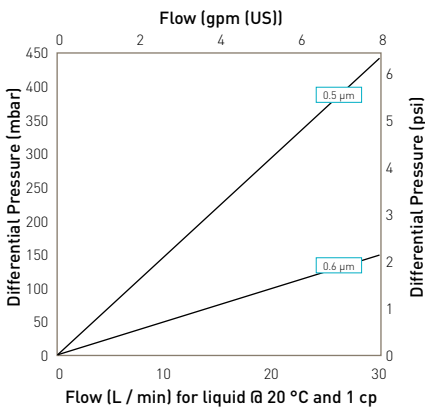

For K size for a given flow rate multiply 10" size differential pressure by 2

10" size (250 mm) filters

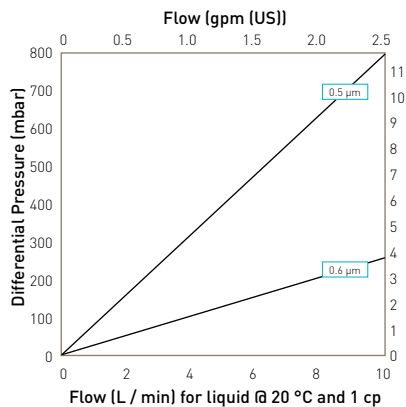

For A size for a given flow rate divide B size differential pressure by 2
For E size for a given flow rate multiply B size differential pressure by 2

B size (125 mm) filters

Specifications

Materials of Construction

- Filtration Media: Glass Microfibre / Polypropylene
- Upstream Support: Polypropylene
- Downstream Support: Polypropylene
- Inner Support Core: Polypropylene
- Outer Protection Cage: Polypropylene
- End Caps: Polypropylene
- End Cap Insert (if applicable): 316L Stainless Steel*
- *Not available in B & L endcap variants
- Standard o-rings/gaskets: Silicone / EPDM
- Capsule Body: Polypropylene
- Capsule Vent Seals: Silicone

Food and Biological Safety

Materials conform to the relevant requirements of 21CFR Part 177, EC1935 / 2004 and current USP Plastics Class VI - 121 °C and ISO10993 equivalents.

Recommended Operating Conditions

Up to 70 °C (158 °F) continuous operating temperature and higher short-term temperatures during CIP to the following limits:

Temperature		Max. Forward dP	
°C	°F	(bar)	(psi)
20	68	5.0	72.5
40	104	4.0	58.0
60	140	3.0	43.5
80	176	2.0	29.0
90	194	1.0	14.5
>100 (steam)	>212 (steam)	0.3	4.0

Capsules may be operated up to a temperature of 40 °C (104 °F) at line pressures up to 5.0 barg (72.51 psig) for liquids.

Effective Filtration Area (EFA)

10" (250 mm) Up to 0.37 m² (3.9 ft²)

Cleaning and Sterilization

PREPOR GP cartridges can be repeatedly steam sterilized in situ or autoclaved at up to 121 °C (249.8 °F). They can be sanitized with hot water at up to 90 °C (194 °F) and are compatible with a wide range of chemicals. Capsules can be repeatedly autoclaved up to 130 °C (266 °F).

Retention Characteristics

The retention characteristics of PREPOR GP have been determined by a combination of controlled laboratory tests and in-use monitoring for a number of organisms. Bacterial challenge testing is carried out to methods specified in ASTM F838-05.

Organism	Approx. Cell Size (µm)*	Typical Titre Reduction			
		0.5	0.6	1.0	1.5
<i>Serratia marcescens</i>	0.5 - 0.8 x 0.9 - 2.0	10 ⁴	10 ³	-	-
<i>Stenococcus oenos</i>	0.5 - 0.7 x 0.7 - 1.2	10 ⁴	10 ³	-	-
<i>Escherichia coli</i>	1.1 - 1.5 x 2.0 - 6.0	10 ⁴	10 ³	-	-
<i>Saccharomyces cerevisiae</i>	1.0 (spherical buds)	10 ⁷	10 ⁶	10 ⁴	10 ³

Ordering Information

Cartridges

ZCGP [] - [] - [] - [] - []

Code Length (Nominal)	Code Micron	Code Endcap (10")	Code Variant	Code O-rings
B* 2.5" (65 mm)	.50 0.5 µm	B* dh DOE	E Electronics	E EPDM
A* 5" (125 mm)	.60 0.6 µm	C BF / 226 Bayonet	S* Steam Sterilizable	P PTFE Encapsulated Silicone
K 5" (125 mm)	.80 0.8 µm	D Fin / 222		S* Silicone
1 10" (250 mm)	1.0 1.0 µm	E Flat Top / 222		V Viton
2 20" (500 mm)	1.5 1.5 µm	G Recess / 222		
3 30" (750 mm)		H UF Retrofit		
4 40" (1000 mm)		J SOE (no o-ring)		
		L* dh DOE		
		N Internal 213		
		R BF / 222 Bayonet		
		Code Endcap (Demi)		
		SK Retrofit		
		T TRUESEAL		
		Y Demi Stub		
		Z Demi A & B Std		

* Supplied in packs of 3.

* For detailed operational procedures and advice on cleaning and sterilization, please contact the Technical Support Group through your usual Parker domnick hunter contact.

* Silicone o-ring supplied as standard without having to specify the 'S' code.

* EPDM gaskets supplied as standard

Capsules

ZEGP [] - [] - [] - [] - [] - [] - [] - [] - []

Code Length (Nominal)	Code Micron	Code Inlet Connection	Code Outlet Connection	Code Variant	Code Grade	Code Pack N°	Code Accessory
E 4.4" (113 mm)	.50 0.5 µm	T 1" Tri-Clamp	T 1" Tri-Clamp	P Pharmaceutical	N Non-Sterile	3 Pack of 3	FB Filling Bell
B 5.5" (140 mm)	.60 0.6 µm	N 1/2" NPT Male	N 1/2" NPT Male				
A 7.9" (200 mm)	.80 0.8 µm	H 1/2" Hosebarb	H 1/2" Hosebarb				
	1.0 1.0 µm	G Stepped Hosebarb	G Stepped Hosebarb				
	1.5 1.5 µm	M 1/4" NPT Male	M 1/4" NPT Male				
		V 3/8" NPT Female	V 3/8" NPT Female				

G & H connections only

* Approx. values as in Holt, J.G., Krieg, N.R., Sneath, P.H.A., Staley, J.T., Williams, S.T., 1994. *Bergey's Manual of Determinative Bacteriology*, Ninth Edition, Williams & Wilkins.

PROSEP FILTER SYSTEMS LTD
Unit G19
River Bank Way
Lowfields Business Park
Elland
West Yorkshire
HX5 9DN

Tel: 01422 377367

Fax: 01422 377369

Email: enquiries@prosep.co.uk

www.prosep.co.uk

Map and Directions to Prosep Filters Limited



Leave M62 at Junction 24.

At roundabout adjacent to Cedar Court Hotel take 2nd exit onto dual carriageway (A629), signposted Halifax.

Take 1st exit slip road.

At roundabout at end of sliproad, take 3rd exit off.

This is the entrance to Lowfields Business Park.

Proceed straight over 1st roundabout.

At next roundabout take 2nd exit onto River Bank Way - Prosep Filters can be found on the left after the S-bend.

[Link to Google Maps](#)